

Texas Tech University

Australian Consumer Panel Line Measuring

- 1. Use Starrett Electronic Digital Caliper.
- 2. Turn Caliper on by pressing **Zero On/Off** button.
- 3. Close Caliper completely and press **Zero On/Off** to calibrate the caliper (make sure reading is in mm).
- 4. To measure lines place inside of stationary prong on the inside of the line labeled 0.
- 5. Slide movable prong until the inside of the prong is in the middle of the line drawn by the consumer.
- 6. (Make sure you measure where the consumer's line intersects the scale).
- 7. Record measurement to the side of the line and initial.
- 8. Each line must be measured by two separate people.
- 9. Make sure caliper stays calibrated by sliding caliper closed completely and making sure reading is 0.00mm.
- 10. When measuring is complete turn off caliper by holding the **Zero On/Off** button until caliper shuts off.
- 11. Place caliper back in case and return it.